



Routine Checks for Extrel Mass Specs

Air Conditioner

- Filter is clean and free of debris
- Analyzer enclosure is between 25°C to 30 °C

Electronics

- All green lights on the 5221

Pumps

- Turbo pump – indicator light is solid green
- Turbo pump – is cool to the touch
- Roughing pump – oil is between level indicators
- Roughing pump – oil is clear/colorless, free of particulates

Sample Entry

- Valve filters are clean or have been replaced within the last 6 months
- Flow switch is clean
- Vacuum pressure is between 1 and 3 x10⁻⁶ Torr

Software

- Questor5 Event Log is free of failures or alarms
- Questor5 Backup databases are less than 75% full



Need a more in-depth assessment?

Extrel's Technical Support team is available to set up a free analyzer performance consultation.

Have questions or need assistance?
Contact us today!

Extrel CMS, LLC

575 Epsilon Drive, Suite 2
Pittsburgh, PA 15238

+1 (412) 967 5754
support@extrel.com
Extrel.com